

Application Data: Adhesives & Sealants Industry

Computrac® MAX® 2000XL Moisture/Solids Analyzer

MATERIAL	RESULT	COMPUTRAC®	STATISTICS	OVEN
Adhesive Resin	% Solids	71.949	Mean	72.087
		0.059	Standard Deviation	0.397
		4 minutes	Test Time	2 hours
Automotive Sealant, Rubber	% Moisture	0.226	Mean	0.228
		0.021	Standard Deviation	0.006
Cold Seal Adhesive	% Moisture	0.544	Mean	0.520
		0.004	Standard Deviation	To constant weight
Glue	% Solids	57.489	Mean	57.586
		0.116	Standard Deviation	0.082
		6 minutes	Test Time	To constant weight
Organic Solvent Base Adhesive	% Solids	33.255	Mean	33.258
		0.158	Standard Deviation	0.158



ARIZONA INSTRUMENT LLC

3375 N. Delaware Street, Chandler, AZ 85225 • (602) 470-1414 • (800) 528-7411 • www.azic.com

Exclusive Manufacturer of Computrac® Moisture Analyzers and Jerome® Toxic Gas Analyzers