

# APPLICATION DATA – FOODS

## Computrac® MAX® 5000

MATERIAL	RESULT	MAX 5000	STATISTIC	OVEN
<b>Beans, BBQ Baked</b> Report #8049	% Moisture	64.0797 0.366 23-25 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	64.3248 0.895 4 hours
<b>Beans, Navy</b> Report #8046	% Moisture	70.9173 0.468 23-25 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	71.6305 0.488 4 hours
<b>Chicken, Deboned</b> Report #7969	% Moisture	70.9796 0.509 8-10 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	70.5185 0.842 18 hours
<b>Chicken, Ground (Whole)</b> Report #7967	% Moisture	72.7452 0.960 10-13 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	72.2509 0.392 18 hours
<b>Egg, Spray Dried</b> Report #7795	% Moisture	4.7493 0.057 2 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	4.7625 0.033 5 hours
<b>Turkey Liver</b> Report #7971	% Moisture	76.6310 0.099 10-13 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	76.6913 0.049 18 hours

MATERIAL	RESULT	MAX 5000	STATISTIC	FURNACE
<b>Beans, Navy</b> Report #8075	% Ash	4.1407 0.067 9-10 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	4.1545 0.028 4 hours
<b>Beer</b> Report #8166	% Ash	0.1104 0.002 21-22 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	0.1015 0.001 3-4 hours
<b>Chicken, Deboned</b> Report #7970	% Ash	6.6759 0.455 16-18 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	6.2324 1.249 2 hours
<b>Chicken, Ground (Whole)</b> Report #7968	% Ash	10.8694 1.233 13-15 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	10.4258 1.039 2 hours
<b>Corn Starch</b> Report # 7786	% Ash	0.1631 0.005 37-38 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	0.1228 0.009 3 hours
<b>Egg, Spray Dried</b> Report #7796	% Ash	4.6997 0.119 31-33 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	4.6932 0.213 4 hours
<b>Turkey Liver</b> Report #7972	% Ash	4.6629 0.219 36-43 minutes	<b>Mean</b> <b>Standard Deviation</b> <b>Test Time</b>	4.6203 0.566 2 hours



**ARIZONA INSTRUMENT LLC**

1912 W. 4<sup>th</sup> Street, Tempe, AZ 85281 • (602) 470-1414 • (800) 528-7411 • www.azic.com

Exclusive Manufacturer of Computrac® Moisture Analyzers and Jerome® Toxic Gas Analyzers