

APPLICATION DATA – PAINT & COATINGS

Computrac® MAX® 5000

MATERIAL	RESULT	MAX 5000	STATISTIC	OVEN
Paint, Water Based Report #6068	% Solids	78.5575 0.101 2 minutes	Mean Standard Deviation Test Time	78.069 0.063 60 minutes
Paint, Solvent Based Report #6067	% Solids	76.1842 0.165 2 minutes	Mean Standard Deviation Test Time	76.8319 0.054 60 minutes
Paint, Stain Primer Sealer Report #7479	% Solids	66.3982 0.422 3 minutes	Mean Standard Deviation Test Time	66.3246 0.554 60 minutes
Paint, Semi Gloss House Report #7473	% Solids	50.9462 0.096 13-14 minutes	Mean Standard Deviation Test Time	51.1347 0.203 60 minutes

MATERIAL	RESULT	MAX 5000	STATISTIC	FURNACE
Paint, Water Based Report #6069	% Ash	76.7448 0.047 5-6 minutes	Mean Standard Deviation Test Time	76.5550 0.101 60 minutes
Paint, Solvent Based Report #6070	% Ash	73.1227 0.085 9-10 minutes	Mean Standard Deviation Test Time	76.3974 0.317 60 minutes
Paint, Stain Primer Sealer Report #7480	% Ash	73.9436 0.096 31-39 minutes	Mean Standard Deviation Test Time	73.4435 0.160 60 minutes
Paint, Semi Gloss House Report #7474	% Ash	51.1795 0.022 10 minutes	Mean Standard Deviation Test Time	50.7653 0.130 60 minutes



ARIZONA INSTRUMENT LLC

1912 W. 4th Street, Tempe, AZ 85281 • (602) 470-1414 • (800) 528-7411 • www.azic.com

Exclusive Manufacturer of Computrac® Moisture Analyzers and Jerome® Toxic Gas Analyzers