

Pharmaceuticals Application Data

Computrac® MAX 2000XL

Material		Computrac	Oven
Emulsion	Mean	47.532% solids	47.487%
	S.D.	0.045%	0.106%
	Test Time	4 minutes	18 hours
Gel Caps (2-piece)	Mean	16.63% moisture	16.62%
	S.D.	0.02%	0.05
	Test Time	8 minutes	17 hours
Granulation	Mean	2.51% moisture	2.53%
	S.D.	0.070%	0.03%
	Test Time	6 minutes	5 hours

Material		Computrac	Titration
Cough Drops	Mean	0.844% moisture	0.814%
	S.D.	0.12%	0.08%
	Test Time	5 minutes	
Dextrose	Mean	8.942% moisture	8.7494%
	S.D.	0.294%	0.163
	Test Time	3 minutes	
Sodium Bicarbonate	Mean	0.01% moisture	0.012%
	S.D.	0.001%	0.002%
	Test Time	3 minutes	
Vancomycin	Mean	11.266% moisture	11.23%
	S.D.	0.021%	0.297%
	Test Time	5 minutes	



ARIZONA INSTRUMENT LLC

1912 W. 4th Street, Tempe, AZ 85281 • (602) 470-1414 • (800) 528-7411 • www.azic.com

Exclusive Manufacturer of Computrac® Moisture Analyzers and Jerome® Toxic Gas Analyzers